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Instruction for Use TPP Centrifuge Tube

The TPP centrifuge tubes are suitable for preparing, mixing, centrifuging and storing suspensions (cells, bacteria, precipitates, etc.) as well as solid and liquid reagents and samples.

The centrifuge tube is intended for single use only.

Note:

Follow national regulations when handling biological material and chemical substances, use appropriate protective clothing

Take note of the regulations of aseptic working during the work process.

Instructions:

- *Check the packaging material and the product for intactness. Use only faultless material.*
- *Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.*
- *Open the package in a sterile environment and remove a complete system for use*
- *Remove the cap and fill according to your laboratory routine. Please refer to the maximal filling quantity (see technical data).*
- *Close the screw cap firm close for gas and liquid tight seal without damaging the sealing lip in the cap. Make sure that the contents cannot escape.*
- *Centrifuging: The use of appropriate rotors and/or centrifuge adapters is recommended. Please follow the safety instructions of the centrifuge manufacturer.*
- *The mechanical load capacity of the centrifuge tubes is influenced by*
 - *Shape and material*
 - *Accuracy of fit in the centrifuge adapter*

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Advice and Recommendations:

- *Centrifuging at (too) high speed can deform the tube.*
- *The graduation serves as a reference for filling quantity and is not an absolute value.*
- *Alcohol 90 % with mechanical stress (rubbing) might dissolve or blur the white color of the print.*

Storage in Minus Temperatures:

- *The product should only be used for its intended use. TPP does not assume any product warranty in case of damage when used otherwise.*
- *For storage in temperatures below °C transfer the samples into TPP cryo tubes. When storing the polypropylene (PP) centrifuge tubes in the minus temperature range, TPP recommends that the product be tested beforehand.*
- *Polystyrene (PS) tubes are not suitable for storage in minus temperatures.*

Important:

- *The material of the tubes and caps is not intended for temperatures minus °C. For storage below °C transfer the samples into TPP cryotubes.*
- *Temperatures below 0 °C result in reduced elasticity for polypropylene. Do not stress frozen tubes mechanically such as dropping or hitting.*

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Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch.