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Instruction for Use TPP Tissue Culture Dish

TPP tissue culture dishes are for the cell/tissue cultivation and growth in manual use. The circumferential gripping ring allows a secure grip of the unit (bowl + lid). The growth surface - but not the side wall of the dish - is opto-mechanically activated for optimal cell adhesion and growth.

The tissue culture dish is for single use.

Note:

Follow national regulations when handling biological material, use appropriate protective clothing

Take note of the regulations of aseptic working during the work process.

Instructions:

- Check the packaging material and the product for intactness. Use only faultless material.
- Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.
- Open the packaging in a sterile environment and remove a complete system (lid with base) for use. The circumferential grip ring allows a secure grip of the cell culture dish and thus avoids unintentional lifting of the lid (risk of contamination).
- Remove the lid and fill the dish with the medium and inoculum according to your laboratory routine.
- To achieve an uniform cell growth, the following aspects must be observed:
 - Resuspend the cells well with the medium to avoid cell lumps.
 - Avoid foaming during resuspension and seeding to avoid holes in the cell layer.
 - Low medium volume leads to meniscus formation and thus to uneven growth of

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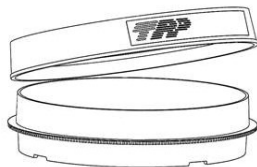
Fig 1, 45 ° abrir

the cells (cell clumps at the edge). Please refer to the technical data for the optimum filling volume.

- *The spacer cams on the inside of the lid ensure a constant and optimal gas exchange with low evaporation.*
- *Tip: For sowing, medium change, sub cultivation and other activities the lid should be opened only ~45° to avoid contamination (Figure 1)*

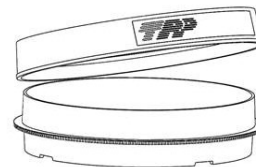
Fig. 1 45° open

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- *Air vents in the bottom rim ensure optimum heat distribution in the incubator even when several dishes are stacked on top of each other.*
- *For identification use the labelling fields.*
- *As an orientation aid, for example during microscopy, the numbers in the clock system 3, 6, 9, 12 are located in the bottom of the dish (Figure 2).*
- *Before use, store at room temperature, protect from UV light*
- *Test the multitude of influencing factors under routine conditions beforehand.*

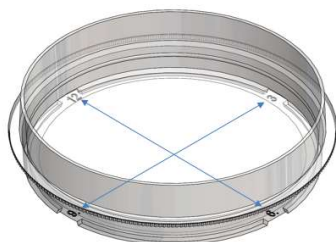
Fig 2, números

Fig. 2, Numbers

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Advice and Recommendations:

- *The stacking rim provides a safe stability of several TPP cell culture dishes stacked on top of each other.*
- *For a homogenous cultivation of the cells, levelling of the incubator and continuous supply of medium is essential.*
- *Avoid and/or eliminate vibration sources near the incubator to ensure uniform cell growth.*
- *Centrifugation of the dish is not recommended.*

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Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch.