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Instructions for Use TPP Syringe Filter

TPP Syringe Filters 22 and 45 are designed for use in filtration of aqueous solutions such as cell culture media and biological fluids of microorganisms, particles and undissolved precipitates larger than 0.22 µm or 0.45 µm - depending on the filter. They can be used bidirectional: Forwards (from the syringe into the container) or backwards (from the container into the syringe), but only in one direction per filter unit and application.

The syringe filter is intended for single use.

Note:

Follow national regulations when handling biological material, use appropriate protective clothing

Take note of the regulations of aseptic working during the work process.

Instructions:

- Check the packaging material and the product for intactness. Use only faultless material.
- Check the expiry date (EXP) on the label and packaging. Use only products with valid EXP.

Forward Filtration:

- Take a syringe and remove the syringe plunger.
- Under sterile conditions, remove the cover paper from the blister packaging of the syringe filter.
- Place the syringe on the Luer-Lock or cone connector of the syringe filter, grasp it while placing it on the blister.

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Reverse Filtration:

- *Remove a syringe from the blister pack.*
- *First draw some air into the syringe to avoid a possible backflow of the liquid, then draw the liquid to be filtered from the container into the syringe.*
- *Remove adhering liquid residues from the syringe tip.*
- *Under sterile conditions, remove the cover paper from the blister packs of the syringe filter.*
- *Place the syringe on the Luer-Lock or cone connector of the syringe filter, grasping it while placing it on the blister.*
- *Filter the liquid by evenly pressing the syringe plunger into a suitable sterile container.*

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Fig1

Remarks and Recommendations:

- *The diameter of the TPP syringe filter fits optimally on a standard 50 ml centrifuge tube. (Fig. 1)*

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Technical Data:

Materials

Membrane Hydrophilic polyether sulfate (PES),
type "TPP fast flow"
Housing PET

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Chemical Compatibility:

Compatible with most aqueous solutions, please consult the relevant documentation on www.tpp.ch.

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Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch.