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### TPP-Dreifach-Verpackung 3-B SAL 10<sup>-6</sup>

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## Instruction for Use

### TPP Triple Bagged 3-B SAL 10<sup>-6</sup>

*The 3-B package is for areas where sterile product transfer is essential. The high Sterility Assurance Level (SAL) of 10<sup>-6</sup> of the products ensures maximum sterility.*

*3-B products are for single use only.*

#### Note:

*Follow the national regulations for handling biological materials and wear the appropriate protective clothing.*

*During the work process, be sure to observe the rules of aseptic technique.*

#### Instructions:

- *Inspect packaging and product for integrity. Use only materials in perfect condition.*
- *Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.*
- *Each product is triple-wrapped in three easy peel-open bags.*
- *Each bag is marked with a security number:  
3 – outer bag,  
2 – middle bag,  
1 – inner bag.*
- *Bag 1 features product data: description, lot number, and EXP.*

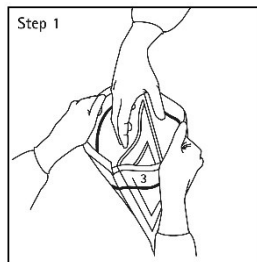
#### Delivery time:

*Please note that this product may not be in stock but produced to order; a delay in delivery may occur.*

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
**Step 1 / Bag 3,  
Entrance Area**

- Open bag 3 in the non-sterile environment.

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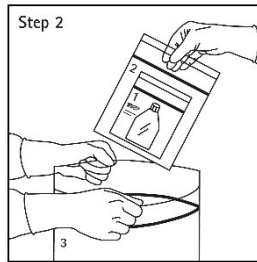
**Step 2 / Bag 2, Transfer**

- Pull out bag 2.
- Transfer to a cleanroom look.

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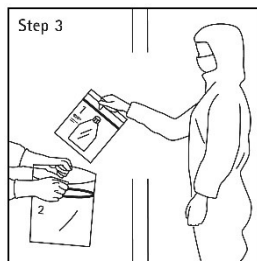**Step 3 / Bag 1, Cleanroom Look**

- Open bag 2 in the cleanroom look or sterile environment, and pull out bag 1.

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**Step 4 / Bag 1, Clean Room**

- Remove product from bag 1 under the biological safety cabinet.

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- For safety, the bag sizes are to prevent damage to the inner bag during opening, thus preventing contamination of the product.

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**Additional:**

- SAL-value:  $10^{-6}$  for high sterilization requirements.
- Full TPP product line available worldwide.

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Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from [www.tpp.ch](http://www.tpp.ch).