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Instructions for Use TPP Tissue Culture Flasks with Peel-off Foil

TPP tissue culture flasks with peel-off foil are for cell culture with the necessary sterile handling techniques, storage and microscopy in manual use and to meet the need for a large one-time access from the above to the cell culture. For optimal cell adhesion and growth, only the bottom of the tissue culture flask is optomechanically activated. The angled neck reduces the risk of medium contacting the inside of the screw cap by spillage and allows excellent access for cell scrapers and serological pipettes.

TPP cell culture flasks with peel-off foil are only available with filter screw caps, if a VENT screw cap is used, a possible pressure build-up of 0.03 bar or higher could push open the peel-off foil.

The TPP tissue culture flask with peel-off foil is for single use only.

Note:

Follow the national regulations for handling biological materials and wear the appropriate protective clothing.

During the work process, be sure to observe the rules of aseptic technique.

Instructions:

- Inspect packaging and product for integrity. Use only materials in perfect condition.
- Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.
- Open the package in a sterile environment and remove a complete system for use.

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Remove the peel-off foil:

- *With the peel-off foil removed, cells can be harvested from the top, i.e., the large opening.*
- *Place your thumb and middle finger on the inscription fields on either side of the of the T-flask neck. Use your index finger to press the flask onto the surface by placing the fingertip underneath the foil flap, thus stabilizing the flask. Grab the foil flap with the other hand and pull the foil back with a gentle pull.*

Important:

- *It is **not** possible to reclose the foil.*

Advice and Recommendations:

- *For labelling there are bilateral inscription fields.*
- *The volume scale is an indication of the filling volume, not an absolute value.*
- *The stacking rim allows a stable stacking of several flasks on top of each other, also in combination with other TPP flasks of the same size.*
- *For homogeneous cultivation of cells, it is essential to level the incubator and provide a continuous supply of medium.*
- *Avoid and/or eliminate sources of vibration near the incubator to ensure uniform cell growth.*

Note:

- *The barrier limits the growth area to 115 cm².*

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Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch.