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Instruction for Use TPP Triple Bagged 3-B

The 3-B packaging is intended for use in areas where the sterile transfer of product is essential. The 3-B solution meets the increasingly more stringent clean room requirements. With the standard single packaging, contamination adhering to the outside of the bag could be unintentionally carried into cleanroom areas. Pouch cleaning is necessary. The 3-B packaging system eliminates this step. The high «Sterility Assurance Level» (SAL) of 10^{-6} of the products guarantees maximum sterility.

The 3-B products are intended for single use only.

Note:

Follow national regulations when handling biological material, use appropriate protective clothing.

Take note of the regulations of aseptic working during the work process.

Instructions:

- *Check the packaging material and the product for intactness. Use only faultless material.*
- *Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.*
- *Each product is triple-wrapped in three easy peel-open bags offering ease of handling*
- *Each bag is marked with a security number:
3 – outer bag
2 – middle bag
1 – inner bag*
- *Bag 1 contains product data: description, lot number and EXP.*

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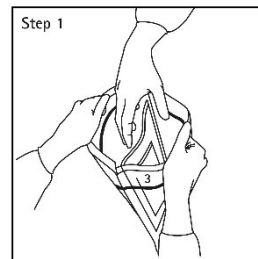
Step 1, bag 3, entrance area

- *Carefully open the outer bag 3 in the non-sterile environment.*
- *Pull out bag 2 by staff within the sterile operating area, ensuring a sterile transfer from one area to the next.*

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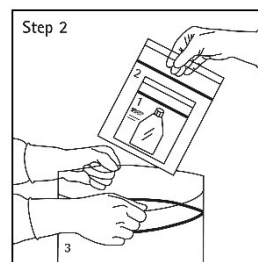
Step 2, bag 2, transfer

- *Pull out bag 2 in sterile environment.*
- *Transfer to sluice*

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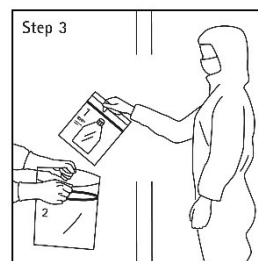
Step 3, bag 1, sluice

- *Open bag 2 in sluice, carefully pull out bag 1 by staff within the sterile operating area, ensuring a sterile transfer through the gowning area to the aseptic zone.*

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Step 4, bag 1, clean room

- *Remove product from bag 1 within the sterile field.*

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- *The safety-open-style peel-open bags are increasing sized, so that during handling of one bag the next inner bag is not accidentally broken rendering it contaminated*

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Additional:

- *SAL-value: 10^{-6} for high sterilization requirements*
- *Entire TPP product range globally available*

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch.

Delivery time: *Note that this product range is not a stock item but manufactured to order, a minimum order quantity may apply.*