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Instruction for Use TPP Tissue Culture Dish

TPP tissue culture dishes are for the cell/tissue cultivation and growth in manual use. The circumferential gripping ring allows a secure grip of the unit (dish + lid). For optimal cell adhesion and growth, only the bottom of the tissue culture dish is opto-mechanically activated.

The tissue culture dish is for single use only.

Note:

Follow the national regulations for handling biological materials and wear the appropriate protective clothing.

During the work process, be sure to observe the rules of aseptic technique.

Instructions:

- *Inspect packaging and product for integrity. Use only materials in perfect condition.*
- *Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.*
- *Open the package in a sterile environment and remove a complete system for use (dish + lid). The circumferential grip ring allows a secure grip of the cell culture dish and thus avoids unintentional lifting of the lid (risk of contamination).*
- *Remove the lid and fill the dish with the medium and inoculate according to your laboratory routine.*
- *To obtain uniform cell growth, the following points must be observed:*
 - *Resuspend the cells well in the medium to avoid cell clumping.*
 - *Avoid foaming during resuspension and seeding to avoid holes in the cell layer.*

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Advice and Recommendations:

- *The stacking rim provides a safe stability of several TPP cell culture dishes stacked on top of each other.*
- *For homogeneous cultivation of cells, it is essential to level the incubator and provide a continuous supply of medium.*
- *Avoid and/or eliminate sources of vibration near the incubator to ensure uniform cell growth.*
- *Centrifugation of the dish is not recommended.*

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Additional:

Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from www.tpp.ch.