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Instructions for Use TPP Tissue Culture Flasks with Re-closable Lid

TPP tissue culture flasks with re-closable are for cell culture with the necessary sterile handling techniques, storage and microscopy in manual use and meet the need for a large one-time access from above to the cell culture. For optimal cell adhesion and growth, only the bottom of the tissue culture flask is opto-mechanically activated. The angled neck reduces the risk of medium contacting the inside of the screw cap by spillage and allows excellent access for cell scrapers and serological pipettes.

TPP tissue culture flasks with peel-off foil are only available with filter screw caps. If a VENT screw cap is used, a possible pressure build-up of 0.03 bar or higher could push open the lid.

The TPP tissue culture flask with re-closable lid is for single use only

Note:

Follow the national regulations for handling biological materials and wear the appropriate protective clothing.

During the work process, be sure to observe the rules of aseptic technique.

Instructions:

- *Inspect packaging and product for integrity. Use only materials in perfect condition.*
- *Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.*
- *Open the package in a sterile environment and remove a complete system for use.*

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Opening Mechanism:

The lid can be opened and closed multiple times and remain gas- and liquid-tight.

- 1 Press here
- 2 Lift out lid

Important: To open the lid, place the flask on a solid surface. To open the re-closable lid, press on the recessed area. Finger pressure on the opening mechanism must be controlled. Excessive pressure may cause breakage.

Advice and Recommendations:

- For labelling there are bilateral inscription fields.
- The volume scale is an indication of the filling volume, not an absolute value.
- The stacking rim allows a stable stacking of several flasks on top of each other, also in combination with other TPP flasks of the same size.
- For homogeneous cultivation of cells, it is essential to level the incubator and provide a continuous supply of medium.
- Avoid and/or eliminate sources of vibration near the incubator to ensure uniform cell growth.
- Always check that the re-closable lid is properly closed before using it for the first time or after opening it.

Note:

- The barrier limits the growth area to 115 cm².

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Additional:

Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from www.tpp.ch.