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## Instruction for Use TPP Cell Scraper and Spatula (Lifter)

*The TPP cell scraper and spatula (or lifter) are used for harvesting adherent cell cultures in various containers such as cell culture dishes, plates and bottles.*

*The angled and pivoting blade simplifies the harvesting procedure and allows for complete cell harvesting with minimal cell damage.*

*The cell scraper and cell lifter is for single use only.*

### Note:

*Follow the national regulations for handling biological materials and wear the appropriate protective clothing.*

*During the work process, be sure to observe the rules of aseptic technique.*

### Instructions:

- *Inspect packaging and product for integrity. Use only materials in perfect condition.*
- *Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.*
- *Open the package in a sterile environment and remove a complete system for use*

### Fields of Application:

- *The cell scrapers are available in 3 sizes to suit your needs (see selection guide).*
- *In addition to traditional cell culture applications, the sterile, individually packaged scrapers and spatulas (lifter) can be used for a variety of other applications.*

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#### Additional:

*Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from [www.tpp.ch](http://www.tpp.ch).*