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## Instructions for Use

### TPP Tissue Culture Flat Tube 10 cm<sup>2</sup> TPP Tissue Culture Tube 20 cm<sup>2</sup>

TPP tissue culture tubes are for the cell/tissue cultivation and growth in manual use. For optimal cell adhesion and growth, only the bottom of the tissue culture dish is opto-mechanically activated.

TPP tissue culture flat tube 10 cm<sup>2</sup> are available with filter screw cap, TPP tissue culture tubes 20 cm<sup>2</sup> are available with VENT screw caps.

The tissue culture tubes are for single use.

#### Note:

Follow the national regulations for handling biological materials and wear the appropriate protective clothing.

During the work process, be sure to observe the rules of aseptic technique.

### Tissue Culture Flat Tube 10 cm<sup>2</sup>, # 91243

#### Instruction:

- Inspect packaging and product for integrity. Use only materials in perfect condition.
- Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.
- Open the package in a sterile environment and remove a complete system for use.
- Cultivate and harvest cells according to your common laboratory routine.
- To obtain uniform cell growth, the following points must be observed:
  - Resuspend the cells well in the medium to avoid cell clumping.

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#### **Filter Screw Cap:**

- *Close the filter screw cap with ventilation holes. Continuous gas exchange occurs through the integrated 0.22 µm hydrophobic membrane. If the membrane becomes wet, gas exchange is temporarily reduced.*
- **Centrifugation:** *The use of appropriate rotors or centrifuge adapters is recommended. Follow the centrifuge manufacturer's safety instructions.*
- *The mechanical strength of the tubes during centrifugation is influenced by*
  - *Shape and material*
  - *Accuracy of fit in the centrifuge adapter*
  - *Temperature, centrifugation time, relative centrifugal force (RCF)*
  - *Chemical and physical properties of the centrifuged material*
  - *Rotor type: fixed-angle or swing-out rotor*
- **Microscopy:** *with inverse microscopes, the flattened top minimizes light refraction.*
- *For identification, use the lateral inscription field. Test the many influencing factors in advance under routine conditions.*
- *Store at room temperature before use, protected from UV light.*

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## Tissue Culture Tube 20 cm<sup>2</sup>, # 91106

### Instruction:

- *Inspect packaging and product for integrity. Use only materials in perfect condition.*
- *Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.*
- *Open the package in a sterile environment and remove a complete system for use.*
- *Cultivate and harvest cells according to your common laboratory routine.*
  
- *To obtain uniform cell growth, the following points must be observed:*
  - *Resuspend the cells well in the medium to avoid cell clumping.*
  - *Avoid foaming during resuspension and seeding to avoid holes in the cell layer*
  - *Low volume of medium leads to meniscus formation and thus to uneven cell growth (cell clumps at the edge). Please refer to the technical data for the optimal filling volume.*

### VENT Screw Cap:

- *For gas exchange, screw the VENT screw cap tight until you hear / feel it clicks into the venting position.*
- *To interrupt the gas exchange, turn the cap a 1/4 turn clockwise over the venting position. The tube is now tightly close there is no gas exchange.*

*For a transfer of the tube within in the laboratory, TPP recommends to close the VENT screw cap tightly.*

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#### **Advice and Recommendations:**

- The volume scale is an indication of the filling volume, not an absolute value.
- For homogeneous cultivation of cells, it is essential to level the incubator and provide a continuous supply of medium.
- Avoid and/or eliminate sources of vibration near the incubator to ensure uniform cell growth.
- PS tubes are not suitable for storage at minus temperatures.

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**Additional:**

Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from [www.tpp.ch](http://www.tpp.ch).