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Instruction for Use TPP Serological Pipettes

Serological pipettes are design for the dispensing of aqueous liquids such as cell culture media or biological liquids.

The color-coding on pipette and dispenser enables easy recognition.

The negative graduation and the patented reservoir provide more working volume than conventional serological pipettes.

The specially designed dispenser with perforated peel-off tab serves as a dispenser for easy collection directly at the laboratory bench.

Note:

Follow national regulations when handling biological material and chemical substances, use appropriate protective clothing

Take note of the regulations of aseptic working during the work process.

Do not use the mouth to pipette; always employ the pipetting-aid to fill and to dispense.

Instructions:

- Check the packaging material and the product for intactness. Use only faultless material.
- Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.
- Open the package in a sterile environment and remove a pipette for use
- Tear open the paper/plastic wrapping to expose the pipet's mouthpiece.
- Place the mouthpiece into the pipetting-aid, keeping the pipette inside its wrapping to avoid contamination. If the pipet does not fit snugly, replace the silicone pipet holder within the pipet aid nozzle.
- With the pipette in position, hold the aid and then remove the wrapping.
- Fill the pipette to the required volume.

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- *Do not overfill the pipette! Avoid liquid coming into contact with the cotton plug or getting into the pipetting aid.*
- *Accuracy is best assured when the pipette is held vertically in front of the eye and the volume measured at the base of the meniscus.*
- *Please consider parallax errors.*
- *Position-the pipette into the target vessel and allow the liquid to run in slowly (on the vessel wall in the case of centrifuge tubes).*
- *Test the multitude of influencing factors under routine conditions beforehand.*
- *Before use, store at room temperature and protect from UV light.*

Advice and Recommendations:

- *If pipetting in a sterile environment and the pipet touches any non-sterile surface, dispose of the pipet and replace it with a sterile one.*
- *Risk of contamination in case of multiple use and/or solution change of the pipette: Always replace the pipette when changing solutions or cell suspensions to avoid cross-contamination.*
- *Always use a pipetting aid (e.g.: Turbo-Fix # 94700, Peleus ball, Aspirette) and never "mouth pipette". For sterile work in cell culture, electronic pipetting aids are the means of choice. All pipetting aids have a sterile filter in the device helps to prevent contamination.*
- *Never place the pipetting aid on the workbench with a filled and/or used pipette. Liquid could flow back into the pipetting aid, contaminate and/or damaging it.*
- *Avoid the formation of aerosols during sterile work with biological agents. Let the pipette drain slowly. Never blow out the last drop out of a mechanical pipette; otherwise, the liquid film at the tip will burst. (Aerosol formation)*
- *Non-sterile solutions can be blown out to the last drop.*
- *Always hold the pipette with the tip down and avoid any contact with your hand.*
- *The serological pipette is not suitable for highly viscous liquids such as glycerol, PEG.*

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- *No liability for being complete, correct and/or up-to-date*
- *No TPP guarantee applies to goods used other than the registered and intended use*

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Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch.

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