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### Tubo para cultura de células 10 cm<sup>2</sup>

### Tubo para cultura de células 20 cm<sup>2</sup>

#### TPP

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### TPP Tissue Culture Flat Tube 10 cm<sup>2</sup>

### TPP Tissue Culture Tube 20 cm<sup>2</sup>

TPP tissue culture tubes are intended for use for the cell/tissue cultivation and growth, in manual use. For an optimal cell adhesion and proliferation, the growth area is opto-mechanically activated.

TPP tissue culture flat tube 10 cm<sup>2</sup> are available with filter screw cap, TPP tissue culture tubes 20 cm<sup>2</sup> are available with VENT screw caps.

The tissue culture tubes are for single use.

#### Note:

Follow national regulations when handling biological material, use appropriate protective clothing

Take note of the regulations of aseptic working during the work process.

### Tissue Culture Flat Tube 10 cm<sup>2</sup>

#### # 91243

#### Instruction:

- Check the packaging material and the product for intactness. Use only faultless material.
- Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.
- Open the packaging in a sterile environment and remove a complete system for use.
- Cultivate and harvest the cells according to the usual laboratory routine.
- In order to achieve uniform growth on the surface, the following aspects must be considered:
  - Resuspend the cells to avoid cell clumps. (Uneven growth)
  - Avoid foaming during resuspension creating holes in the cell lawn.

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#### **Filter Screw Cap:**

- *For the continuous gas exchange, close the filter screw cap with integrated hydrophobic PTFE filter membrane 0.22 µm tightly. If the membrane gets wet, the gas exchange reduces for a short period.*
- *For identification, use the lateral inscription field.*
- **Centrifuging:** *The use of suitable rotors or centrifuge adapters is recommended. Follow the safety instructions of the centrifuge manufacturer.*  
*The mechanical strength of the tubes during centrifugation is influenced by*
  - *Shape and material*
  - *Accuracy of fit in the centrifuge adapter*
  - *Temperature, centrifugation time, g-number*
  - *Chemical and physical properties of the centrifuged material*
  - *Rotor type: fixed-angle or swing-out rotor*
- **Microscope** *with inverse microscopes, the flattened top minimizes light refraction.*
- *Test the multitude of influencing factors under routine conditions beforehand.*
- *Storage before use at room temperature, protected from UV light.*

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## **Tissue Culture Tube 20 cm<sup>2</sup> # 91106**

- *Check the packaging material and the product for intactness. Use only faultless material.*
- *Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.*
- *Open the packaging in a sterile environment and remove a complete system for use.*
- *Cultivate and harvest the cells according to the usual laboratory routine.*
- *In order to achieve uniform growth on the surface, the following aspects must be considered:*
  - *Resuspend the cells to avoid cell clumps. (Uneven growth)*
  - *Low medium volume leads to the formation of minuscule cells, which leads to an uneven growth of the cells. (many cells on the edge, few in the middle). For optimum filling volume refer to the technical data*

### **VENT Screw Cap:**

- *For gas exchange, screw the VENT screw cap tight until you hear/feel it clicks into the venting position.*
- *To interrupt the gas exchange, turn the cap a 1/4 turn clockwise over the venting position. The tube is now tightly close there is no gas exchange.*

*For a transfer of the tube within in the laboratory, TPP recommends to close the VENT screw cap tightly.*

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#### **Advice and Recommendations:**

- *The graduation serves as a reference for filling quantity and is not an absolute value.*
- *For a homogenous cultivation of the cells, levelling of the incubator and continuous supply of medium is essential.*
- *Avoid and/or eliminate vibration sources near the incubator to ensure uniform cell growth.*
- *PS tubes are not suitable for storage at minus temperatures*

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**Additional:**

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage [www.tpp.ch](http://www.tpp.ch).